SafetyCaps

 **Installationsanleitung**

 **Installation Manual**

Distributed by:

Lab Unlimited
CARL STUART GROUP

Tallaght Business Park
Whitestown, Dublin 24,
Ireland
D24 RFK3

Tel: (01) 4523432
Fax: (01) 4523967
E-mail: info@labunlimited.com
Web: www.labunlimited.com

Quatro House, Frimley Road,
Camberley,
United Kingdom
GU16 7ER

Tel: 08452 30 40 30
Fax: 08452 30 50 30
E-mail: info@labunlimited.co.uk
Web: www.labunlimited.co.uk

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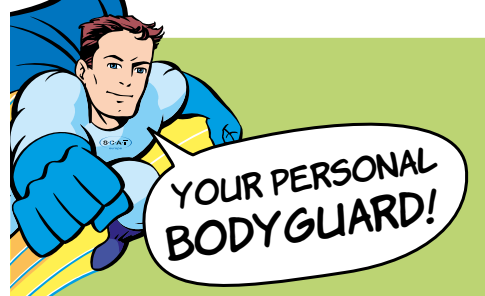
PDF

YOUR PERSONAL BODYGUARD!

- ✓ **Health protection and workspace safety:** Hazardous vapours are reduced by 70%.
- ✓ **Save money:** Prevent evaporation and leakage of expensive solvents.
- ✓ **Improve chromatographic results:** Prevent retention time shifts due to concentration changes, caused by evaporation.
- ✓ **Easy solvent bottle exchange:** No twisting tubes during bottle exchange, thanks to the freely rotating screw cap.

EN 

How to install your SafetyCap:

- 1 Screw the **air valve** into a free port of the SafetyCap.
- 2 Guide the **solvent tube** through the **fitting** first, and then through the **SafetyCap**.
- 3 Reconnect the **solvent filter** to the solvent tube (at the open end).
- 4 Screw the **SafetyCap** on top of your solvent bottle.
- 5 Adjust the **length** of the **solvent tube** in the solvent bottle.
- 6 **Fix** the tube by tightening the **fitting**.

Air valves secure your work environment and are permanently exposed to vapours, dust particles and pollutants. Renew used air valves after 6 months. For an optimal degree of safety.

PDF